

ADICTYE™ TISSUE COLLECTION KIT

INSTRUCTIONS FOR USE



www.adicyte.com

All components of the AdiCyte Adipose Tissue Collection are sterile. Please maintain sterile technique during collection steps to ensure a sterile sample for processing. Extra syringe caps and Luer-Lok™ adapters are included for your convenience or use with other systems.

Collecting from a STERILE Canister/filter: (Common for 60cc syringe kits)

1. Wipe Luer-Lok port on canister/filter with alcohol or betadine wipe.
2. Withdraw syringe barrel to completely fill syringe.
3. Remove syringe from Luer-Lok and seal with sterile syringe cap.
4. Place bar code label on syringe.
5. Place capped syringe in plastic bag. Remove excess air in bag and seal.
6. Place bagged syringe in kit and repeat process for additional syringes.

Collecting with a Cannula: (Common for smaller 10 cc syringe kits)

1. Connect sterile cannula to syringe Luer-Lok.
2. Withdraw syringe barrel to completely fill syringe.
3. Remove syringe from cannula and seal with sterile syringe cap.
4. Place bar code label on syringe.
5. Place capped syringe in plastic bag. Remove excess air in bag and seal.
6. Place bagged syringe in kit and repeat process for additional syringe.

*Tissue may also be transferred from larger to smaller syringes using sterile technique.

AFTER COLLECTION

1. Complete the **Data Collection Sheet**. Place a bar code label in the bottom right corner of the document.
2. Close kit, seal shipping container and **Complete Sender Information on pre-printed Shipping Label**.
3. Adhere shipping label to container.

COURIER PICK-UP INSTRUCTIONS (Samples should ideally arrive within 24 hours of collection)

1. Call FedEx™ Express Priority Overnight at **1-800-463-3339**.
2. Please arrange for first available, same-day pick up. If same-day pick up is not available, the shipment may be dropped off at any FedEx location.
3. ACCOUNT NUMBER: **319913351**
4. Lab Address (already pre-printed)

Dr. David Harris: AdiCyte Cell Bank
University of Arizona Health Sciences Center 6129
1501 North Campbell Ave. Tucson, AZ 85724
Lab Phone: 520-626-1610 Emergency Only 1-888-262-0387

RETURN ALL UNUSED COMPONENTS AND BAR CODES